

pH-Fix 6.0–10.0

Test strips for measuring pH values in the range of 6.0–10.0

 Please refer to instructions for use

REF 92122

Pack content:

100 test strips
1 reference color scale pH 6.0–10.0

Instructions for use:

1. Dip test strip for 1 s with all test fields in the sample solution.
2. Shake off excess sample solution.
3. Compare with color scale and read off the corresponding pH-value.



Note:

If sample solution is weakly buffered leave test strip in solution until color no longer changes.

Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container in a dry place (storage temperature 4–30 °C).

Disposal:

Used test strips may be disposed of as domestic waste.

 Please read instructions for use!

REF Item number

 Manufacturer